NMRCheck : This program checks the existence of new Bruker NMR files on mora every 10 minutes. If something is found, it will be copied to your predefined target folder (e.g. I:\nmr)!

written by: Jan P. Unsleber

## 1) Installation:

- a) Create new folder: e.g. on drive I:\nmr (alternative a local folder or a folder on your group disk)
- b) Download nmrcheck.zip from <u>\\mora\nmr\mnova</u> to C:\Users\<your account>
- c) Click with right mouse button on the zip-file and choose "Extract all" or "Alle extrahieren"

## 2) Start:

- a) Change to the installation directory (e.g. C:\Users\<your account>\nmrcheck)
- b) Click with right mouse button on nmrcheck link and choose "Pin to task bar" or "An Taskleiste anheften"
- c) Click the nmrcheck icon on the task bar
- 3) Choose your settings Hint: By clicking on the buttons, you can toggle between active (white) or inactive (red).

# Main settings:

- a) Choose the right institute, e.g.
- b) Input your **personal** usertag, e.g.

# Advanced settings:

Open Folder

De eluvre Cresseture	If you enable this option, you've to create and input the path of your target folder,
Backup Spectra	where all you're NMR data should be copied to, e.g. I:\nmr

Open in MNova If you enable this option, the copied data is loaded into MNova.

OC

ocf

Path to MNova, **NWZ:** N:\Software\Chemie\MNova\MNova.lnk

local: C:\Program Files (x86)\Mestrelab Research S.L\MestReNova\MNova.exe

If you enable this option, the new NMR data is shown by the explorer.

Display pop-up If you enable this option, a pop-up appears, after new data was found.

#### 4) Start checking/Check now [Enter]

C NMRCheck v	3.0.0				Local tar	get fold	er I:\nmr:		
#	NMR	CHECK	# <u>^</u> 	4					x
   v3.0.0 (:	11.05.2015)	by Jan P. Un	sleber		▶ nmr ▶	<b>▼</b> 4 <sub>7</sub>	nmr durchsuchen		٩
Last check Options rea Checking [] NEW: \\morr NEW: \\morr Checking []	was at [Fri Jun ad from "config. Mon Jul 6 16:31 a\nmr\300er\Jun1 A\nmr\300er\Jun1 Mon Jul 6 17:05	19 11:53:31 2015]. ini". :13 2015] 9-2015\590 9-2015\591 :27 2015]	#		Neuer Ordner	N;	ame ocf_p_70_v_24_33 ocf_p_70_v_24_33	€_Jun19-2015_591	0
	Ma	iin settings:							
Insitute:	ICPC	OC							
Usertag:	ocf								
NMR folder:	\\mora\nmr\300er								
	\\mora\nmr\400er								
	\\mora\nmr\dpx300								
	Adva	nced settings:			Choose a	target	older with v	write perm	issions
Backup spectra	l arget folder:	l:\nmr							_
Open in MNova	Path to MNova:	N:\Software\Chemie\MNova\MI	Nova.Ink		Don't for	get to ir	iput the righ	nt extensio	n
Op	oen Folder	Display pop-up							
	Start checkir	ig/Check now [Enter]			You can o new data be carrie program	click on a instant d out ev is runni	this button a ly! (Otherw rery 15 minu ng)	at any time ise the che ıtes, as lon	e to get ck will g as the