Date: 23.02.2021

Authorized user: T. Grottke, Work area: Cleanroom,

**Physics Lab** 

Workplace: CeNTech I, 00.17

## Safety Instruction Critical Point Dryer Leica EM CPD 300

WWU
Münster Nanofabrication
Facility (MNF)
Center for NanoTechnology
(CeNTech)
Heisenbergstraße 11
48149 Münster

Range of Application	
Working with Critical Point Dryer Leica EM CPD 300	
Hazard for People and the environment	
<ul> <li>Contact with electricity or potential for electric shock</li> <li>Exposure to CO2 gas</li> <li>Exposure to pressure vessel</li> </ul>	
Safety Measures and Rules of Behavior	
The machine is to be used only by instructed persons  General:  It is forbidden to dry any biological samples or any samples that might contaminate this machine  Only the provided ethanol is to be used as exchange medium  Only the provided holder and filler are to be used  The lid may only be removed when the pressure gauge reads 1bar  Turn off instrument, and be sure NOT to tighten nuts on lid, so that the next user knows upon starting work that the dryer is not under pressure  The dryer requires the use of liquid CO2 cylinder with a purity of 5.0. Only trained staff is permitted to move and install these cylinders  Personal protective equipment:  Safety goggles  Lab coat  Nitrile gloves  Enclosed footwear  Storage:	
The provided ethanol is to be stored at a well-ventilated place	
Behavior in a Dangerous Situation	
<ul> <li>Enter the danger zone only with adequate protective clothing</li> <li>Get people out of the danger zone</li> <li>Immediately leave the danger zone</li> <li>Inform your superior</li> <li>In case of fire, every fire-extinguishing medium can be used</li> </ul>	
Action in Accident, First Aid	
<ul> <li>Treat smaller injuries</li> <li>Inform first aiders</li> <li>Note accident in the accident book</li> <li>In case of sever injuries, inform an emergency doctor</li> </ul>	
Inform your superior     Disposal	
Do not dispose of the waste in the drain and use only the provided container.	
To not dispose of the waste in the drain and use only the provided container.	Dücksoital